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- 주조용 알루미늄 (Cast Aluminum Alloy)

Al 3 20% Si 50X BF

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Table 1. Cutting Parameters

NEW TechCut 5x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Silicon Carbide
Wheel Bond	Resin
Size	8"
Wheel RPM	3500
Feed Rate	12.7 mm/min

Table 2. Hot Mounting Parameters

TechPress3™	
Mold Size	1.25" (32 mm)
Powder	Black glass-filled epoxy powder
Pre-Heat	Off
Heat Temp/Time	160°C / 9 min
Cool Time	6 min
Pressure	4.5 bar

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Table 3. MetPrep3™/PH-3™ Gridding/Polishing Procedure

Step	1	2	3	4	5	6	
Consumables	Abrasive Size (P-Grading)	180 Grit (P-180)	320 Grit (P-400)	600 Grit (P-1200)	6 μm	1 μm	0.04 μm
	Type	Silicon Carbide	Silicon Carbide	Silicon Carbide	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Abrasive Disc	Abrasive Disc	Abrasive Disc	Glycol Suspension	Glycol Suspension	Suspension
	Polishing Cloth	-	-	-	Gold Label	DiaMat	Final A
	Coolant	Water	Water	Water	GreenLube	GreenLube	Water**
Settings	Platen Speed(RPM) / Direction	300 Comp	300 Comp	300 Comp	150 Comp	150 Comp	150 Comp
	Sample Speed (RPM)	150	150	150	150	150	150
	Force (lbf/N) +	4/18	4/18	4/18	5/22	5/22	4/18
	Time (min)	1*	1	1	3	2	1

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

Table 4. AD-5™ Dispensing Parameters

Step	4&5	6
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	2.0 sec
Water Flush	Off	On

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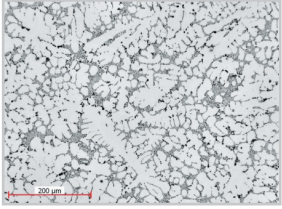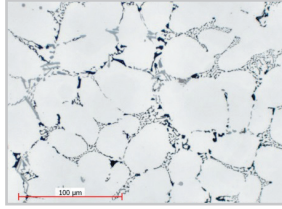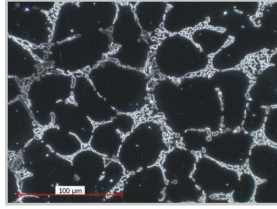
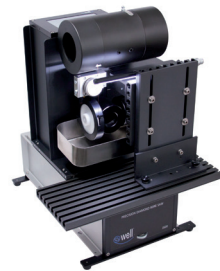
NEW TechCut 5x™ Precision High Speed Saw

WELL 3500 Premium Diamond
Wire Cutting System

TechPress3™ Hydraulic-Pneumatic Mounting
Press / 1 1/4" Mold Assembly

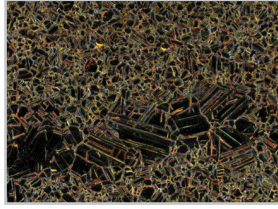
MetPrep3™ / PH-3™ Grinding/ Polishing
System, AD-5™ Automatic Fluid Dispenser

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Bronze 500x DF

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Table 1. Cutting Parameters

PowerCut 10x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Silicon Carbide
Wheel Bond	Resin
Size	10"

Table 2. Hot Mounting Parameters

TechPress3™	
Mold Size	1.5" (38 mm)
Powder	Black glass-filled epoxy powder
Pre-Heat	Off
Heat Temp/Time	165°C / 10 min
Cool Time	7 min
Pressure	5.0 bar

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Table 3. MetPrep3™/PH-3™ Gridding/Polishing Procedure

Step	1	2	3	4	5	6	
Consumables	Abrasive Size (P-Grading)	180 Grit (P-180)	320 Grit (P-400)	600 Grit (P-1200)	6 μm	1 μm	0.04 μm
	Type	Silicon Carbide	Silicon Carbide	Silicon Carbide	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Abrasive Disc	Abrasive Disc	Abrasive Disc	Glycol Suspension	Glycol Suspension	Suspension
	Polishing Cloth	-	-	-	Gold Label	Vel-Cloth	Final A
	Coolant	Water	Water	Water	RedLube	RedLube	Water**
Settings	Platen Speed(RPM) / Direction	300 Comp	300 Comp	300 Comp	150 Comp	150 Comp	150 Comp
	Sample Speed (RPM)	150	150	150	150	150	150
	Force (lbf/N) +	4/18	4/18	4/18	6/27	4/18	4/18
	Time (min)	1*	1	1	2	2	2

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

Table 4. AD-5™ Dispensing Parameters

Step	4&5	6
Type	Diamond Suspension & RedLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	2.0 sec
Water Flush	Off	On

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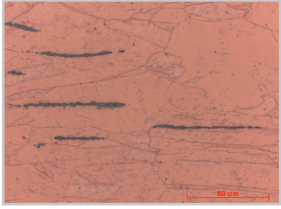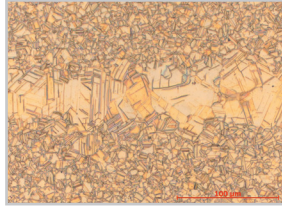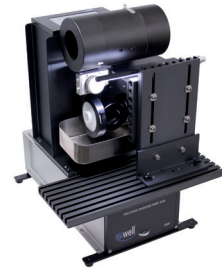
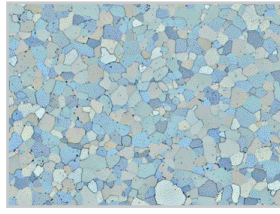
Titanium 100x pol

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Table 1. Cutting Parameters

PowerCut 10x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Silicon Carbide
Wheel Bond	Resin
Size	8"
Wheel RPM	3800
Feed Rate	12.7 mm/min

Table 2. Hot Mounting Parameters

TechPress3™	
Mold Size	1.25" (32 mm)
Powder	Black glass-filled epoxy powder
Pre-Heat	Off
Heat Temp/Time	160°C / 9 min
Cool Time	6 min
Pressure	4.5 bar

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Table 3. MetPrep4™ / PH-6™ Grinding/Polishing Procedure

Step	1	2	3	4	5	
Consumables	Abrasive Size (P-Grading)	180 Grit (P-180)	320 Grit (P-400)	9 μm	3 μm	0.04 μm
	Type	Silicon Carbide	Silicon Carbide	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Abrasive Disc	Abrasive Disc	Glycol Suspension	Glycol Suspension	Suspension
	Polishing Cloth	-	-	PLAN-B	DiaMat	Final A
	Coolant	Water	Water	GreenLube	GreenLube	Water**
Settings	Platen Speed(RPM) / Direction	300 Comp	300 Comp	150 Comp	150 Comp	150 Comp
	Sample Speed (RPM)	150	150	150	150	150
	Force (lbf/N) +	4/18	4/18	5/22	5/22	4/18
	Time (min)	1*	1	3	3	1

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

Table 4. AD-5™ Dispensing Parameters

Step	3&4	5
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	2.0 sec
Water Flush	Off	On

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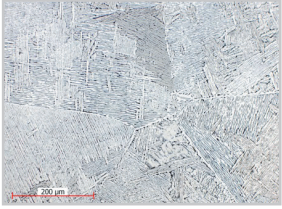
8 Ti-6Al-4V, etched, 200x, Brightfield

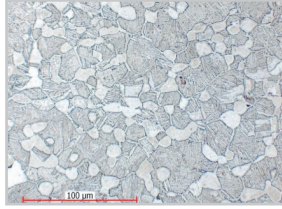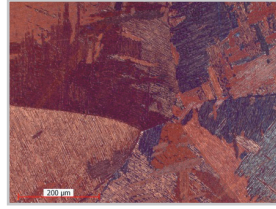Alpha-Beta Titanium, etched, 500x, Brightfield

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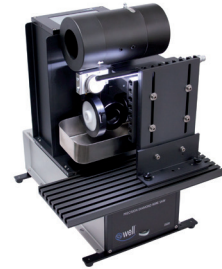
NEW TechCut 5x™ Precision High Speed Saw

WELL 3500 Premium Diamond Wire Cutting System

TechPress3™ Hydraulic-Pneumatic Mounting Press / 1 1/4" Mold Assembly

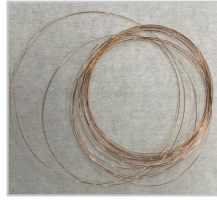
MepPrep4™ / PH-3™ Grinding/ Polishing System, AD-5™ Automatic Fluid Dispenser

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Table 1. How to EpoxyBond 110

Mixed	Part A : Part B = 10 : 1 (By Volume)
Curing Schedule	5 minutes at 150°C 10 minutes at 120°C 30 minutes at 100°C

Table 2. Cold Mounting Parameters

EpoxySet	
Mold Size	1.25" (32 mm)
Mix Ratio (weight) Resin : Hardener	100:12
Cure Time	8Hours (25°C)

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
Table 3. MetPrep4™ / PH-6™ Grinding/Polishing Procedure

Step	Grinding				Polishing	
	1	2	3	4	5	6
Abrasive	320Grit (P-400)	600Grit (P-1200)	800Grit (P-2400)	1200Grit (P-4000)	3 μm	0.04 μm
Type	SiC	SiC	SiC	SiC	Diamond	Colloidal Silica
Carrier	Abrasive Disc	Abrasive Disc	Abrasive Disc	Abrasive Disc	Suspension	Suspension
Polishing Cloth	-	-	-	-	DiaMat	Chem-POL
Coolant	Water	Water	Water	Water	GreenLube	-
Platen Speed(RPM) / Direction	100/Comp	100/Comp	100/Comp	100/Comp	100/Contra	100/Contra
Sample Speed (RPM)	90	90	90	90	90	90
Force (lbf)	3	3	4	4	4	4
Time (min)	Until Target	Until Target	2'00"	2'30"	2'00"	2'00"


Table 4. AD-5™ Dispensing Parameters

Step	5	6
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	6
Pulse Length	0.5 sec	1 sec
Water Flush	Off	On

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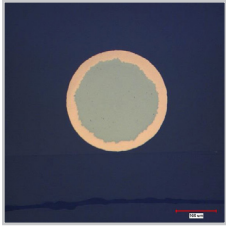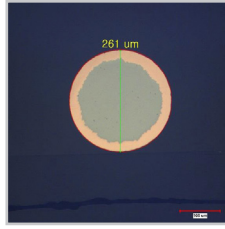
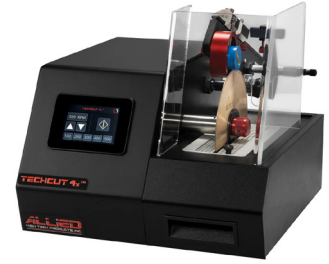
NEW TechCut 4x™ Precision Low Speed Saw

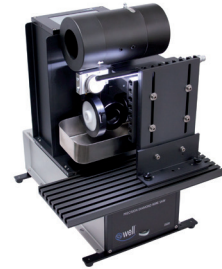WELL 3500 Premium Diamond Wire Cutting System

VacuPrep™ Epoxy Impregnation System

DualPrep3™/PH-4™ Grinding/Polishing System, AD-5™ Automatic Fluid Dispenser

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