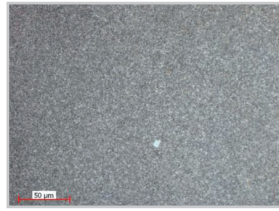
16

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Tungsten carbide, etched, 50x, Brightfield

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Table 1. Cutting Parameters

NEW TechCut 5x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Aluminum Oxide
Wheel Bond	Resin
Size	8"
Wheel RPM	4000
Feed Rate	12.7 mm/min

Table 2. Hot Mounting Parameters

TechPress3™	
Mold Size	1.25" (32 mm)
Powder	Black glass-filled epoxy powder
Pre-Heat	Off
Heat Temp/Time	160°C / 9 min
Cool Time	6 min
Pressure	4.5 bar

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Table 3. MetPrep3™/PH-3™ Gridding/Polishing Procedure

Step	1	2	3	4	5	6	
Consumables	Abrasive Size (P-Grading)	70 µm	15 µm	6 µm	3 µm	1 µm	0.05 µm
	Type	Diamond	Diamond	Diamond	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Metal Bond Dia-Grid	Metal Bond Dia-Grid	Metal Bond Dia-Grid	Glycol Suspension	Glycol Suspension	Suspension
	Polishing Cloth	-	-	-	Gold Label	White Label	Final A
	Coolant	Water	Water	Water	GreenLube	GreenLube	Water**
Settings	Platen Speed(RPM) / Direction	150 Comp	150 Comp	150 Comp	150 Contra	150 Contra	150 Contra
	Sample Speed (RPM)	150	150	150	150	150	150
	Force (lbf/N) *	5/22	5/22	5/22	6/27	6/27	3/13
	Time (min)	2*	2	2	3	2	1

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

Table 4. AD-5™ Dispensing Parameters

Step	4&5	6
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	2.0 sec
Water Flush	Off	On

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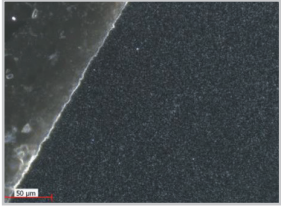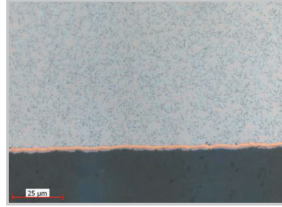Tungsten carbide, as-polished, 50x, Darkfield

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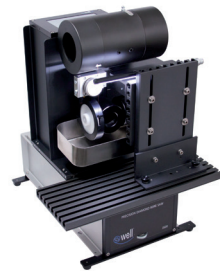
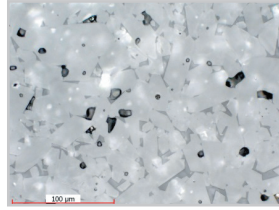
- 알루미나 (Alumina, Al₂O₃)

Aluminum oxide, as-polished, 500x, Brightfield

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Table 1. Cutting Parameters

NEW TechCut 5x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Diamond
Wheel Bond	Resin
Size	7"
Wheel RPM	4000
Feed Rate	12.7 mm/min

Table 2. Cold Mounting Parameters

EpoxySet	
Mold Size	1.25" (32 mm)
Mix Ratio (weight) Resin : Hardener	100 : 12
Cure Time	8 Hours (25°C)

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Table 3. MetPrep3™/PH-3™ Gridding/Polishing Procedure

Step	1	2	3	4	5	6	
Consumables	Abrasive Size (P-Grading)	70 µm	15 µm	9 µm	6 µm	1 µm	0.04 µm
	Type	Diamond	Diamond	Polycrystalline Diamond	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Metal Bonded DiaGrid Disc	Metal Bonded DiaGrid Disc	Suspension	Suspension	Suspension	Suspension
	Polishing Cloth	-	-	PLAN-Cloth	Gold Label	TECH-Cloth	Final A
	Coolant	Water	Water	GreenLube	GreenLube	GreenLube	Water**
Settings	Platen Speed(RPM) / Direction	150 Comp	150 Comp	150 Comp	150 Comp	150 Comp	150 Comp
	Sample Speed (RPM)	150	150	150	150	150	150
	Force (lbf/N) *	5	5	6	6	5	4
	Time (min)	2*	12	4	4	2	1

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

Table 4. AD-5™ Dispensing Parameters

Step	4&5	6
Type	Diamond Suspension & RedLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	2.0 sec
Water Flush	Off	On

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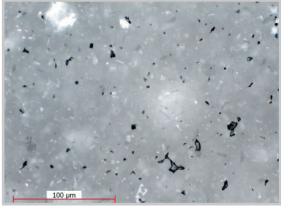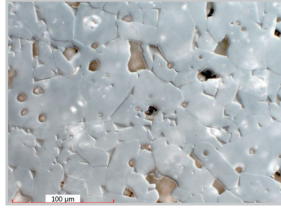Aluminum oxide, as-polished, 500x, Brightfield

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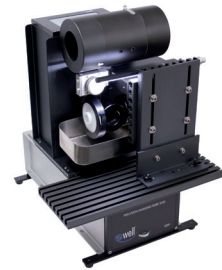
Cement Clinker BF

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Table 1. Cutting Parameters

NEW TechCut 5x™	
Coolant	Lubri-Cut Cutting Fluid
Wheel Type	Diamond
Wheel Bond	Resin
Size	7"
Wheel RPM	3000
Feed Rate	6.35 mm/min

Table 2. Cold Mounting Parameters

EpoxyMount	
Mold Size	1.25" (32 mm)
Mix Ratio (weight) Resin : Hardener	10 : 3
Cure Time	2 Hours (25°C)

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Table 3. MetPrep3™/PH-3™ Grinding/Polishing Procedure

Step	1	2	3	4	5	
Consumables	Abrasive Size (P-Grading)	320 Grit (P-400)	600 Grit (P-1200)	6 μm	1 μm	0.04 μm
	Type	Silicon Carbide	Silicon Carbide	Polycrystalline Diamond	Polycrystalline Diamond	Colloidal Silica
	Carrier	Abrasive Disc	Abrasive Disc	Glycol Suspension	Glycol Suspension	Suspension
	Polishing Cloth	-	-	White Label	White Label	Final A
	Coolant	Water	Water	GreenLube	GreenLube	Water**
Settings	Platen Speed(RPM) / Direction	300 Comp	300 Comp	150 Contra	150 Contra	150 Contra
	Sample Speed (RPM)	150	150	150	150	150
	Force (lbf/N) *	5/22	5/22	5/22	5/22	5/22
	Time (min)	1*	2	8	8	2

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

Table 4. AD-5™ Dispensing Parameters

Step	3&4	5
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	10
Pulse Length	0.5 sec	1.0 sec
Water Flush	Off	On

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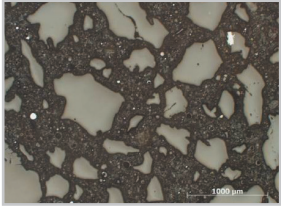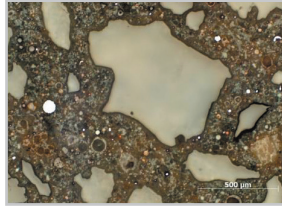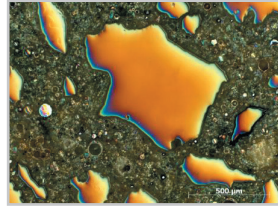
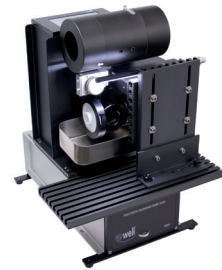
Table 1. Cold Mounting Parameters

EpoxyMount	
Mold Size	1.25" (32 mm)
Mix Ratio (weight) Resin : Hardener	10:3
Cure Time	2Hours (25°C)

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
Table 2. Grinding/Polishing - Auto Polisher

Step	Grinding		Polishing		
	1	2	3	4	5
Abrasive	600Grit (P-1200)	800 Grit (P-2400)	9 μm	3 μm	0.04 μm
Type	Dia-Grid	SiC	Diamond	Diamond	Colloidal Silica
Carrier	Diamond Disc	Abrasive Disc	Suspension	Suspension	Suspension
Polishing Cloth			Gold Label	DiaMat	Chem-Pol
Coolant	Water	Water	GreenLube	GreenLube	-
Platen Speed(RPM) / Direction	100/Comp	100/Comp	100/Contra	100/Contra	100/Contra
Sample Speed (RPM)	90	90	90	90	90
Force (IbF)	4	5	5	5	4
Time (min)	Until Flat	3'00"	5'00"	3'00"	3'00"

Table 3. AD-5™ Dispensing Parameters

Step	3&4	5
Type	Diamond Suspension & GreenLube	Colloidal Silica Suspension
Pulses/Min	6	6
Pulse Length	0.5sec	1sec
Water Flush	Off	On

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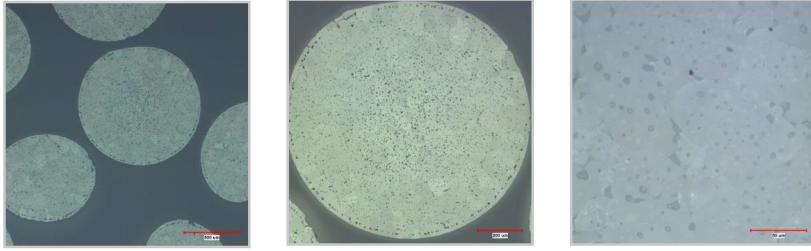
- #145-10005 EpoxyMount Kit
- #197-10005 1¼" 2-Part Mounting Cups
- #50-70805M 600 Grit(P-1200) Dia-Grid Grinding Disc
- #50-10076 800 Grit Silicon Carbide Abrasive Disc
- #90-30030 9 µm Polycrystalline Diamond Suspension
- #90-30020 3 µm Polycrystalline Diamond Suspension
- #180-25000 0.04 µm Colloidal Silica Suspension
- #90-209000 GreenLube™ Polishing Lubricant
- #90-700-210 Gold Label Rigid Back Disc
- #90-700-550 DiaMat Rigid Back Disc
- #180-10750 Chem-Pol Rigid Back Disc
- #148-10000 Micro Organic Soap

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
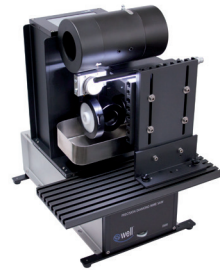
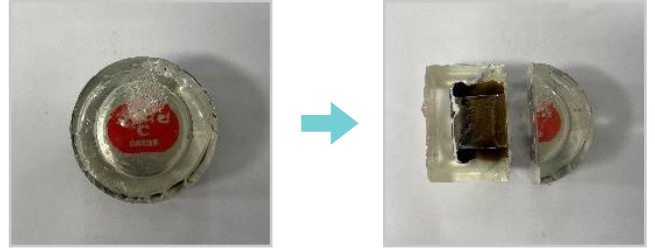
Table 2. AD-5™ Dispensing Parameters

Step	5
Type	Diamond Suspension & GreenLube
Pulses/Min	6
Pulse Length	0.5sec
Water Flush	Off

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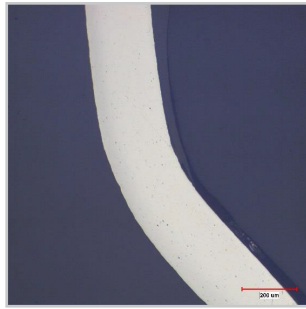
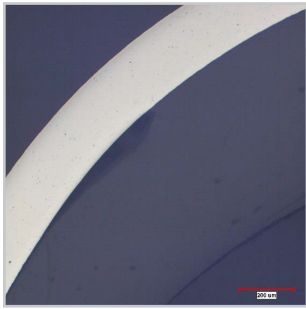
Glass

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Table 1. Cold Mounting Parameters

EpoxyMount	
Mold Size	1.25" (32 mm)
Mix Ratio (weight) Resin : Hardener	100:12
Cure Time	8Hours (25°C)

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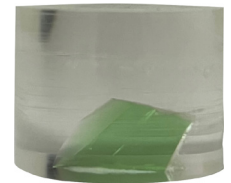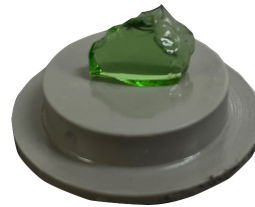Table 2. Grinding/Polishing - Auto Polisher

Step	Grinding		Polishing	
	1	2	3	4
Abrasive	600Grit (P-1200)	800Grit (P-2400)	3 μm	0.5 μm
Type	SiC	SiC	Diamond	Cerium Oxide
Carrier	Abrasive Disc	Abrasive Disc	Suspension	Suspension
Polishing Cloth	-	-	DiaMat	Final A
Coolant	Water	Water	GreenLube	-
Platen Speed(RPM) / Direction	80/Comp	80/Comp	80/Comp	80/Comp
Sample Speed (RPM)	60	60	60	60
Force (lbF)	3	4	3	3
Time (min)	Until Flat	2'30"	2'00"	1'00"


Table 3. AD-5™ Dispensing Parameters

Step	3
Type	Diamond Suspension & GreenLube
Pulses/Min	6
Pulse Length	0.5sec
Water Flush	Off

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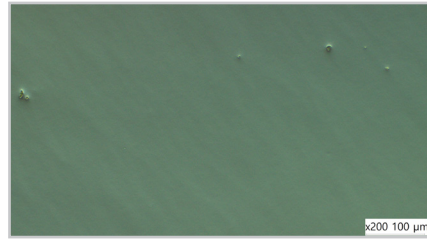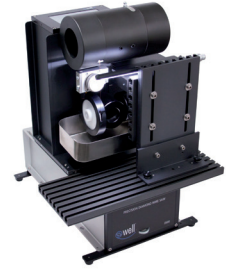
1. 비정형유리
2. Quartz Glass Blocks
3. Soda-Lime Glass Discs
4. Silicate Glass Disc
5. Quartz Glass Fiber Optic Rods

WELL 3500 Premium Diamond Wire Cutting System

NEW TechCut 5x™ Precision High Speed Saw

VacuPrep™ Epoxy Impregnation System

MetPrep3™ / PH-3™ Grinding/ Polishing System, AD-5™ Automatic Fluid Dispenser

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